



LONGGENE OPTIMAL SOFTWARE INSTRUCTIONS FOR USE

In order to use your Real Time PCR devices effectively and smoothly, please follow the instructions below and follow them completely;

1. USE A VIRUS FREE FAT32 FORMAT USB DISK (32GB OR LESS) TO IMPORT PROTOCOLS OR EXPORT YOUR RUN FILES FROM THE UNIT.
2. TO UPLOAD A PROTOCOL TO INSTRUMENT (TXC FILE):
 - a. PREPARE YOUR 'TXC' FILE BY USING A COMPUTER
 - b. COPY IT TO USB DISK AND PLUG IT INTO THE FRONT USB PORT OF THE INSTRUMENT.
 - c. MAIN MENU> ALL PROGRAM> USB DISK> SELECT YOUR PROTOKOL (TXC FILE)
 - d. FILES> COPY> PUBLIC FOLDER (SELECT IT FROM THE DROP DOWN MENU)
 - e. SELECT AND RUN YOUR PROTOCOL BY USING> ALL PROGRAM> PUBLIC> PROTOCOL NAME (TXC FILE)
3. TO DOWNLOAD A RUN FILE FROM THE INSTRUMENT (ZSJ FILE):
 - a. PLUG YOUR USB DRIVE TO FRONT SIDE OF THE INSTRUMENT
 - b. MAIN MENU> QPCR DATA> SELECT YOUR FILE> EXPORT
 - c. PLUG THE USB DRIVE TO YOUR COMPUTER
 - d. SELECT YOUR FILE (ZSJ), COPY IT TO YOUR COMPUTER
 - e. OPEN OPTIMAL SOFTWARE, SELECT OPEN> YOUR FILE> RESULTS.